



BATTERY INSTALLATION AND REPLACEMENT

A. Battery installation position: Right bottom of the freezer body and left outside the electrical box

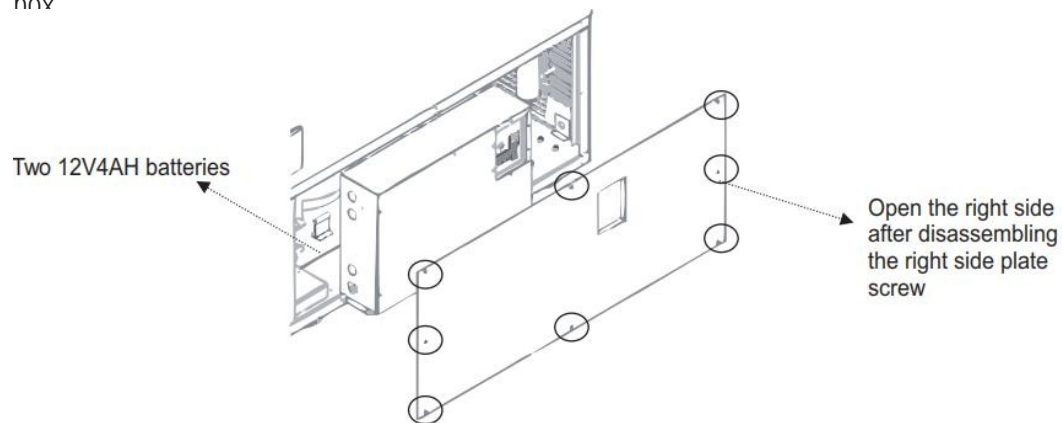

B. Replace the battery:

- Turn off the power switch and pull out the power plug from the outlet (Note: There are high-voltage electrical components in the electrical box, therefore before opening, turn off the power supply, unplug the power plug and turn off the power switch of the freezer, and the electrical box can only be opened by qualified engineers or maintenance personnel).
- Remove the screws on the right-side plate with a screwdriver. Unplug the battery and connect the plug cord. Before unplugging the plug cord, pay attention to the cathode/anode of the battery and the sequence of the connecting line, so as to prevent burning down the control system due to installation of the new battery with cathode/anode errors. Generally, the red wire is connected to the cathode, and the black wire is connected to the anode.

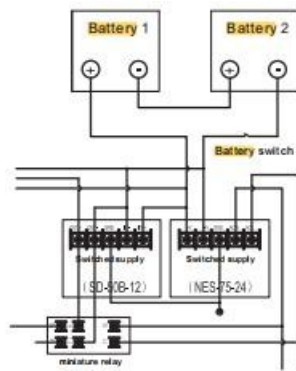

d. Remove the battery fixing clip and take out the battery.

e. The newly replaced batteries shall be two 12V/4AH batteries.

f. Battery disposal:

For a replaced battery that can be recycled, please contact the local battery recycling agency. Note: In order to effectively ensure that the battery replacement of the freezer meets the requirements of the control system and avoid the influence of improper operation on the system during the replacement process, it is recommended to contact the after-sales personnel of Kalstein France for replacement or guidance.

